



Use of 2-5°C cold storage facilities

Humidity is kept at 20-30%.

1671-015: Intended for smaller soil or sediment material and water samples.

1674-013: Intended for larger material and sediment cores.

1. Cooling facilities are for **short term storage of geological samples only**.
2. Keep the room and the entrance **tidy and clean**. Spills on shelves or floors must be removed immediately.
3. **Use plastic boxes for all smaller items**, no single/loose samples, bags or bottles.
4. All **boxes and larger sample items** should be **labeled with date, name and contact info of the owner** using a marker. For external samples, please add the **name and contact info of the responsible employee** to the label. Markers, laminated sheets and extra bags are stored at the entrance.
5. **No wood, paper or cartridge boxes, as well as pieces of paper with information**, they support mold.
6. Use **air-tight containers or wrap your items in material of suitable size** to avoid cross-contamination between samples.
7. Any irregularities regarding access, temperature, humidity etc. must be reported to the lab manager.

Samples store incorrectly are discarded, without further notice. The status of the cooling facilities is checked monthly.